


FLEXCELL®

INTERNATIONAL CORPORATION

JULY 2017 NEWSLETTER

Vol. 4
No. 3

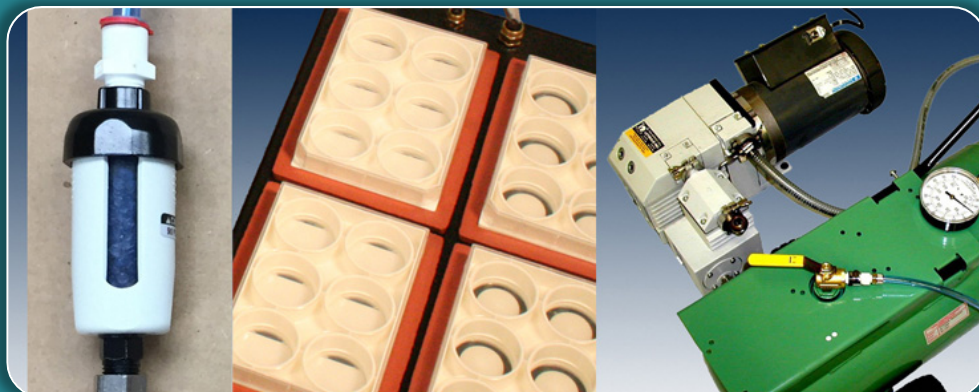

SALE!

5% OFF*

FX-5000™ Tension System
FX-5000™ Compression System
Tissue Train® System

**Sale ends July 31, 2017*

Upcoming Tradeshows

ASCB/EMBO 2017
December 2-6, 2017
Philadelphia, PA

*See website each month for
more details*

www.flexcellint.com

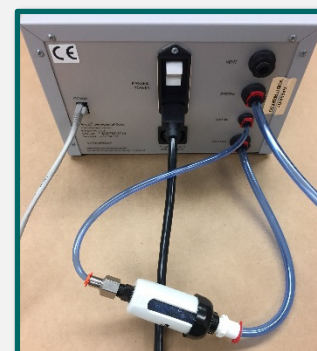


Figure 1. Drying Filter in-line with FX-5000™ Tension System.

Preventative Maintenance

Keep your Flexcell® equipment and components in better working order with regular preventative maintenance. We recommend the following:

- Use the **drying filter** (Fig. 1) whenever the system has been operated for more than one day and will not be operated again for more than one day. This routine maintenance will prolong the life of the valves and reduce expensive repairs.*
- Clean baseplates and gaskets with regular laboratory detergent, rinse with warm tap water followed by distilled or deionized water, and allow to dry.
- Store baseplates flat. Ensure that the entire bottom of baseplate is supported to prevent warping.
- Store gaskets on a flat surface (do not hang) and away from UV light exposure.
- Store culture plates at room temperature or 4 °C in the dark or out of direct light.
- Change the oil on your Leybold Vacuum Pump after the first 100 operating hours and then every 2000-3000 operating hours or after one year. Oil should appear transparent and yellowish in color.
- Flush flow devices and tubing with dH₂O to ensure all culture medium or flow perfusate is removed. Remove pressure on silicone tubing when not in use.
- See user manuals for more information on maintaining Flexcell® equipment.

* Failure to run the drying regimen as recommended can void the warranty.



Medi Analytika

Meet Medi Analytika...

Distributor for Flexcell® International in India... Medi Analytika is a leading instrumentation company in India with over 20 years of experience. Learn more about them at www.medianalytika.com.

Flexcell® International Corp. • 2730 Tucker Street, Suite 200 • Burlington, NC 27215
(919) 732-1591 • (800) 728-3714 • info@flexcellint.com • www.flexcellint.com